


Recommended Use Of The Terra Core®



NOTE: The Terra Core® Sampler is a single use device. It cannot be cleaned and/or reused.

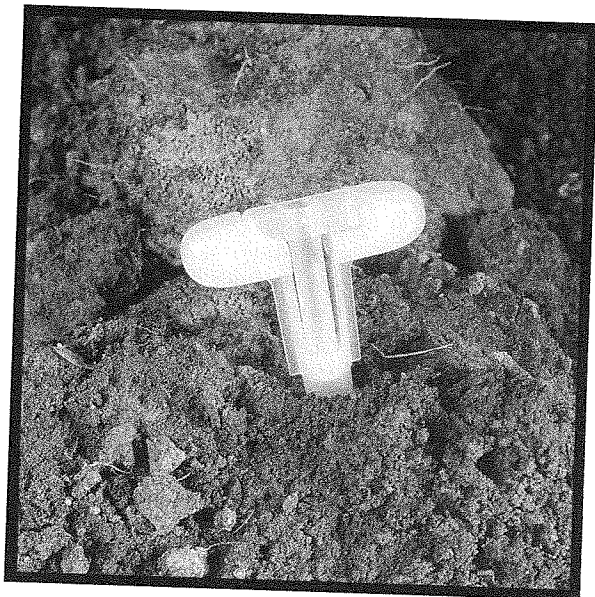


Step 1

Have ready a 40ml glass VOA vial containing the appropriate preservative. With the plunger seated in the handle, push the Terra Core® into freshly exposed soil until the sample chamber is filled. A filled chamber will deliver approximately 5 or 10 grams of soil.

Step 2

Wipe all soil or debris from the outside of the Terra Core® sampler. The soil plug should be flush with the mouth of the sampler. Remove any excess soil that extends beyond the mouth of the sampler.

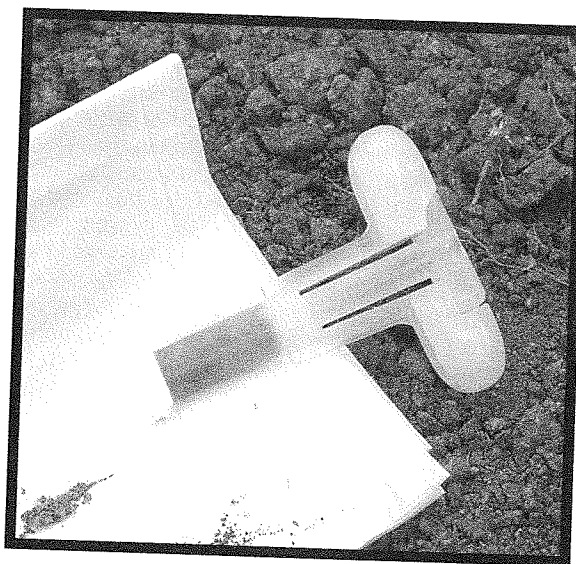

Step 3

Rotate the plunger that was seated in the handle top 90° until it is aligned with the slots in the body. Place the mouth of the sampler into the 40ml VOA vial containing the appropriate preservative and extrude the sample by pushing the plunger down. Quickly place the lid back on the 40ml VOA vial. **Note:** When capping the 40ml VOA vial, be sure to remove any soil or debris from the top and/or threads of the vial.

